




Bedienungsanleitung

Adapterscheibe **GeccoFix (Flex)**

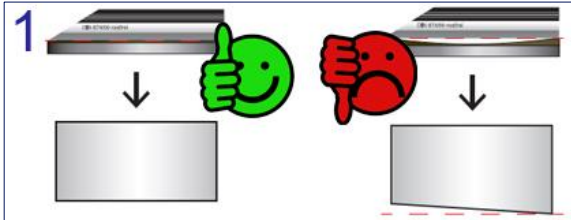


1 Vorbereitung


Stellen Sie sicher, dass die **Arbeitscheibe** Ihrer Maschine (inkl. **Magnetadapter**) **sauber und plan** ist. Gegebenenfalls Planlauf mit geeignetem Messmittel prüfen.

2 **GeccoFix (Flex)** auflegen

GeccoFix (Flex) auf die Arbeitscheibe legen und anschließend die **Schutzfolie der Haftseite** entfernen.

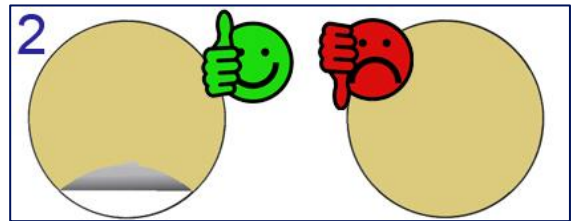

Hinweis: Bei **selbstklebendem Verbrauchsmaterial** wird empfohlen, die **Schutzfolie auf dessen Rückseite nicht zu entfernen**.


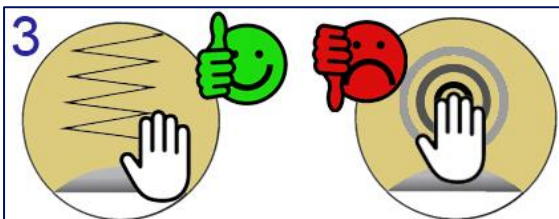


3 Verbrauchsmaterial anbringen

Schleifpapier/Poliertuch/Schleifolie an **einem Punkt anlegen** und anschließend in **Zickzack-Bewegung** über das Material streichen, damit es **blasenfrei** aufliegt.

4 Betrieb

Während des Schleif-/Polierprozesses sollte der **Probenbeweger leicht außerhalb** der **GeccoFix (Flex)** positioniert sein.

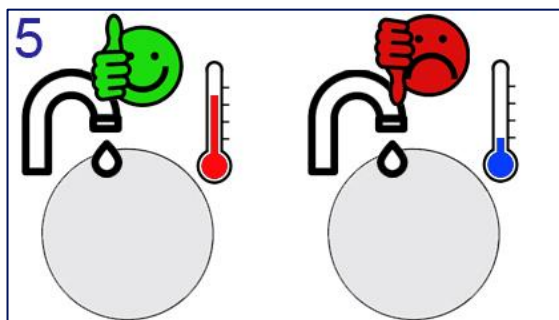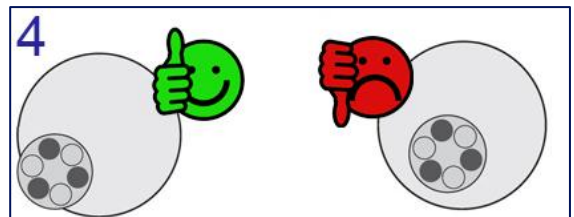

5 Reinigung

GeccoFix (Flex) mit **lauwarmem Wasser** und etwas **Spülmittel** reinigen. Vor erneutem Gebrauch mit **fusselfreiem Tuch** gründlich **trocknen** oder **an der Luft trocknen** lassen.




EXTRA: Auf **GeccoFix (Flex)** können **normales Schleifpapier**, **foliertes Papier** sowie **alle Arten selbstklebender Tücher & Papiere** und **Schleiffolien** verwendet werden.



USER GUIDE


adapter disk **GeccoFix (Flex)**

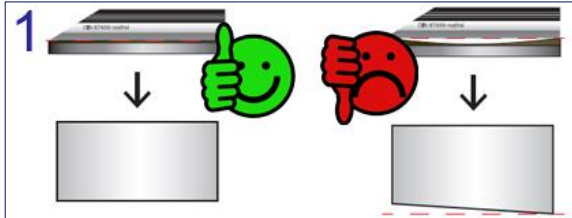


1 Preparation

Make sure the **machine platen** (incl. magnetic adapter) is **clean and flat**. Check flatness with suitable measuring equipment if needed.

2 Place **GeccoFix (Flex)**
Lay **GeccoFix (Flex)** on the platen, then peel off the **liner** from its adhesive side.

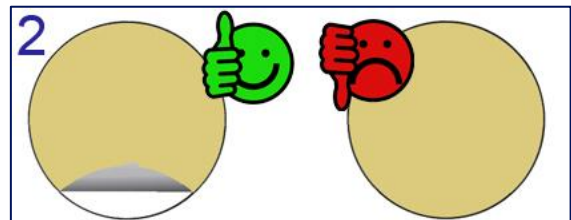



Note: When using **self-adhesive** consumables, it is recommended **not** to remove the liner on their back side.

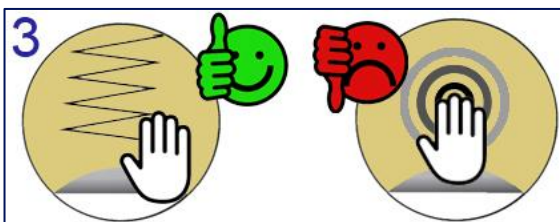

3 Apply the consumable

Position the grinding paper / polishing cloth / grinding foil **at one point** and **press it on in a zigzag motion** to **avoid air bubbles**.

4 Operation
During grinding/polishing process, the **specimen mover** should be positioned **slightly outside** the edge of the **GeccoFix (Flex)**.

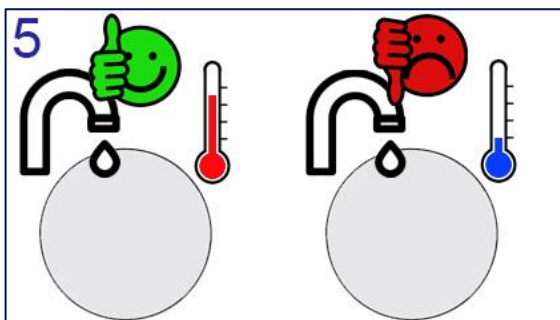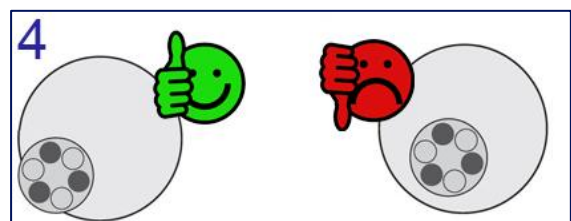

5 Cleaning

Clean **GeccoFix (Flex)** with **lukewarm water and mild dish soap**. Before reuse, **dry thoroughly** with a lint-free cloth or let it air-dry.



EXTRA: **GeccoFix (Flex)** can be used with **standard grinding papers, backed/foiled papers, and all types of self-adhesive cloths & papers** as well as **grinding foils**.